

# Anti-Phospho-Ser<sup>845</sup> GluR1 Antibody



**PhosphoSolutions®**

Antibodies that work™

www.phosphosolutions.com

order@phosphosolutions.com

888-442-7100

**Catalog #:** p1160-845

**Size:** 150 µl

**Cite this Antibody:** PhosphoSolutions Cat# p1160-845, RRID:AB\_2492128

Host	Applications	Species Tested	Species Reactivity*	Molecular Weight
Rabbit	WB 1:1000 IHC 1:1000 (frozen sections)	M, R, turtle	H, NHP	~100 kDa

**Product Description:** Affinity purified rabbit polyclonal antibody.

**Biological Significance:** The ion channels activated by glutamate are typically divided into two classes. Those that are sensitive to N-methyl-D-aspartate (NMDA) are designated NMDA receptors (NMDAR) while those activated by  $\alpha$ -amino-3-hydroxy-5-methyl-4-isoxalone propionic acid (AMPA) are known as AMPA receptors (AMPA). The AMPAR are comprised of four distinct glutamate receptor subunits designated (GluR1-4) and they play key roles in virtually all excitatory neurotransmission in the brain (Keinänen et al., 1990; Hollmann and Heinemann, 1994). The GluR1 subunit is widely expressed throughout the nervous system. Phosphorylation of Ser<sup>845</sup> on GluR1 is thought to be mediated by PKA and phosphorylation of this site increases the conductance of the AMPAR (Roche et al., 1996; Banke et al., 2000). In addition, phosphorylation of this site has been linked to synaptic plasticity as well as learning and memory (Lee et al., 2003; Esteban et al., 2003).

**Antigen:** Phosphopeptide corresponding to amino acid residues surrounding the phospho-Ser<sup>845</sup> of rat GluR1.

**Antibody Specificity:** Specific for endogenous levels of the ~100 kDa GluR1 protein phosphorylated at Ser<sup>845</sup>. Immunolabeling is completely eliminated by treatment with  $\lambda$ -Ptase.


**Purification Method:** Prepared from pooled rabbit serum by affinity purification via sequential chromatography on phospho and non-phosphopeptide affinity columns.

**Quality Control Tests:** Western blots performed on each lot.

**Packaging:** 150 µl in 10 mM HEPES (pH 7.5), 150 mM NaCl, 100 µg BSA per ml and 50% glycerol.

**Storage and Stability:** Shipped on blue ice. Storage at -20°C is recommended, as aliquots may be taken without freeze/thawing due to presence of 50% glycerol. Stable for at least 1 year at -20°C.

**Product Specific References:**

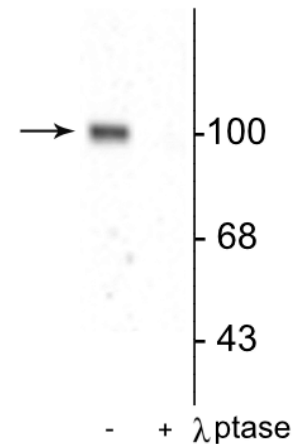

Western blot of rat hippocampal lysate showing specific immunolabeling of the ~100 kDa GluR1 protein phosphorylated at Ser<sup>845</sup> in the first lane (-). Phosphospecificity is shown in the second lane (+) where immunolabeling is completely eliminated by blot treatment with  $\lambda$  phosphatase ( $\lambda$ -Ptase, 1200 units for 30 min).

- Hikida, T., Morita, M., Kuroiwa, M., Macpherson, T., Shuto, T., Sotogaku, N., Niwa, M., Sawa, A. and Nishi, A., 2019. Adolescent psychosocial stress enhances sensitization to cocaine exposure in genetically vulnerable mice. *Neuroscience research*, pii: S0168-0102(19)30021-5.
- Mao, L.M., He, N., Jin, D.Z. and Wang, J.Q., 2018. Regulation of Phosphorylation of AMPA Glutamate Receptors by Muscarinic M4 Receptors in the Striatum In vivo. *Neuroscience*, 375, pp.84-93.
- Suzuki, K., Harada, A., Suzuki, H., Capuani, C., Ugolini, A., Corsi, M. and Kimura, H., 2018. Combined treatment with a selective PDE10A inhibitor TAK-063 and either haloperidol or olanzapine at subeffective doses produces potent antipsychotic-like effects without affecting plasma prolactin levels and cataleptic responses in rodents. *Pharmacology Research & Perspectives*, 6(1).
- Xue, B., Mao, L.M., Jin, D.Z. and Wang, J.Q., 2017. Pharmacological modulation of AMPA receptor phosphorylation by dopamine and muscarinic receptor agents in the rat medial prefrontal cortex. *European journal of pharmacology*. 2017 Dec 11;820:45-52.
- Tronci, E., Napolitano, F., Muñoz, A., Fidalgo, C., Rossi, F., Björklund, A., Usiello, A. and Carta, M., 2017. BDNF over-expression induces striatal serotonin fiber sprouting and increases the susceptibility to l-DOPA-induced dyskinesia in 6-OHDA-lesioned rats. *Nov;297:73-8. Experimental Neurology*.
- Musante, V., Li, L., Kanyo, J., Lam, T.T., Colangelo, C.M., Cheng, S.K., Brody, A.H., Greengard, P., Le Novere, N. and Nairn, A.C., 2017. Reciprocal regulation of ARPP-16 by PKA and MAST3 kinases provides a cAMP-regulated switch in protein phosphatase 2A inhibition. *Elife*, 6, p.e24998.
- Lameth, J., Gervais, A., Colin, C., Lévêque, P., Jay, T.M., Edeline, J.M. and Mallat, M., 2017. Acute Neuroinflammation Promotes Cell Responses to 1800 MHz GSM Electromagnetic Fields in the Rat Cerebral Cortex. *Neurotoxicity Research*, pp.1-16.
- Andrade, E.C., Musante, V., Horiuchi, A., Matsuzaki, H., Brody, A.H., Wu, T., Greengard, P., Taylor, J.R. and Nairn, A.C., 2017. ARPP-16 is a striatal-enriched inhibitor of protein phosphatase 2A regulated by microtubule-associated serine/threonine kinase 3 (Mast 3 kinase). *Journal of Neuroscience*, pp.4559-15.
- Hollis, F., Sevelinges, Y., Grosse, J., Zanoletti, O. and Sandi, C., 2016. Involvement of CRFR1 in the basolateral amygdala in the immediate fear extinction deficit. *eNeuro*, 3(5).
- Xue, B., Fitzgerald, C. A., Jin, D. Z., Mao, L. M., & Wang, J. Q. (2016). Xue, B., Fitzgerald, C.A., Jin, D.Z., Mao, L.M. and Wang, J.Q., 2016. Amphetamine elevates phosphorylation of eukaryotic initiation factor 2 $\alpha$  (eIF2 $\alpha$ ) in the rat forebrain via activating dopamine D1 and D2 receptors. *Brain research*, 1646, pp.459-466.
- Xue, B., Chen, E. C., He, N., Jin, D. Z., Mao, L. M., & Wang, J. Q. (2017). Xue, B., Chen, E.C., He, N., Jin, D.Z., Mao, L.M. and Wang, J.Q., 2017. Integrated regulation of AMPA glutamate receptor phosphorylation in the striatum by dopamine and acetylcholine. *Neuropharmacology*, 112, pp.57-65.
- O'Leary, H., Bernard, P. B., Castano, A. M., & Benke, T. A. (2016). Enhanced long term potentiation and decreased AMPA receptor desensitization in the acute period following a single kainate induced early life seizure. *Neurobiology of disease*, 87, 134-144.
- Tassin, T. C., Benavides, D. R., Plattner, F., Nishi, A., & Bibb, J. A. (2015). Regulation of ERK Kinase by MEK1 Kinase Inhibition in the Brain. *Journal of Biological Chemistry*, Jun 26;290(26):16319-29.
- Mao, L. M., Xue, B., Jin, D. Z., & Wang, J. Q. (2015). Dynamic increases in AMPA receptor phosphorylation in the rat hippocampus in response to amphetamine. *Journal of neurochemistry*. Jun;133(6):795-805.
- Ip, F. C., Fu, W. Y., Cheng, E. Y., Tong, E. P., Lok, K. C., Liang, Y., Ye, W.C. & Ip, N. Y. (2015). Anemoside A3 Enhances Cognition through the Regulation of Synaptic Function and Neuroprotection. *Neuropsychopharmacology*, 40, 1877-1887.
- Suzuki K, Harada A, Shiraishi E, Kimura H. (2015) In Vivo Pharmacological Characterization of TAK-063, a Potent and Selective Phosphodiesterase 10A Inhibitor with Antipsychotic-Like Activity in Rodents. *JPET*, 352:471-479.
- Mango, D., Bonito-Oliva, A., Ledonne, A., Cappellacci, L., Petrelli, R., Nisticò, R., Berretta, N., Fisone, G. and Mercuri, N.B., 2014. Adenosine A1 receptor stimulation reduces D1 receptor-mediated GABAergic transmission from striato-nigral terminals and attenuates l-DOPA-induced dyskinesia in dopamine-denervated mice. *Experimental neurology*, 261, pp.733-743.
- Squillace, M., Doderio, L., Federici, M., Migliarini, S., Errico, F., Napolitano, F., Krashia, P., Di Maio, A., Galbusera, A., Bifone, A. and Scattoni, M.L., 2014. Dysfunctional dopaminergic neurotransmission in asocial BTBR mice. *Translational psychiatry*, 4(8), p.e427.
- Wang, J., Zhang, Y., Xu, H., Zhu, S., Wang, H., He, J., Zhang, H., Guo, H., Kong, J., Huang, Q., & Li, X. M. (2014). Fluoxetine Improves Behavioral Performance by Suppressing the Production of Soluble  $\beta$ -Amyloid in APP/PS1 Mice. *Current Alzheimer Research*, 11(7), 672-680.
- Meyer, D. A., Torres-Altora, M. I., Tan, Z., Tozzi, A., Di Filippo, M., DiNapoli, V., Plattner, F., Kansy, J.W., Benkovic, S.A., Huber, J.D., Miller, D.B., Greengard, P., Calabresi, P., Rosen, C.L., & Bibb, J. A. (2014). Ischemic stroke injury is mediated by aberrant Cdk5. *The Journal of Neuroscience*, 34(24), 8259-8267.
- Mao, L. M., Hastings, J. M., Fibuch, E. E., & Wang, J. Q. (2014). Propofol selectively alters GluA1 AMPA receptor phosphorylation in the hippocampus but not prefrontal cortex in young and aged mice. *European journal of pharmacology*, 738, 237-244.
- Bernard, P. B., Castano, A. M., Bayer, K. U., & Benke, T. A. (2014). Necessary, but not sufficient: insights into the mechanisms of mGluR mediated long-term depression from a rat model of early life seizures. *Neuropharmacology*, 84, 1-12.

- Xue, B., Edwards, M. C., Mao, L. M., Guo, M. L., Jin, D. Z., Fibuch, E. E., & Wang, J. Q. (2014). Rapid and sustained GluA1 S845 phosphorylation in synaptic and extrasynaptic locations in the rat forebrain following amphetamine administration. *Neurochemistry international*, 64, 48-54.
- Shuto, T., Kuroiwa, M., Koga, Y., Kawahara, Y., Sotogaku, N., Toyomasu, K. and Nishi, A., 2013. Acute effects of resveratrol to enhance cocaine-induced dopamine neurotransmission in the striatum. *Neuroscience letters*, 542, pp.107-112.
- Shen Y, Fu WY, Cheng EY, Fu AK, and Ip NY (2013) Melanocortin-4 receptor regulates hippocampal synaptic plasticity through a protein kinase A-dependent mechanism. *J Neurosci*. 33(2):464-72.
- Santini E, Feyder M, Gangarossa G, Bateup HS, Greengard P, Fisone G. (2012) Dopamine- and cAMP-regulated phosphoprotein of 32-kDa (DARPP-32)-dependent activation of extracellular signal-regulated kinase (ERK) and mammalian target of rapamycin complex 1 (mTORC1) signaling in experimental parkinsonism. *J Biol Chem*. 287(33):27806-12.
- Errico, F., Bonito-Oliva, A., Bagetta, V., Vitucci, D., Romano, R., Zianni, E., Napolitano, F., Marinucci, S., Di Luca, M., Calabresi, P. and Fisone, G., 2011. Higher free D-aspartate and N-methyl-D-aspartate levels prevent striatal depotentiation and anticipate L-DOPA-induced dyskinesia. *Experimental neurology*, 232(2), pp.240-250.
- Napolitano, F., Bonito-Oliva, A., Federici, M., Carta, M., Errico, F., Magara, S., Martella, G., Nistico, R., Centonze, D., Pisani, A. and Gu, H.H., 2010. Role of aberrant striatal dopamine D1 receptor/cAMP/protein kinase A/DARPP32 signaling in the paradoxical calming effect of amphetamine. *Journal of Neuroscience*, 30(33), pp.11043-11056.
- Robert M. Sears, Rong-Jian Liu, Nandakumar S. Narayanan, Ruth Sharf, Mark F. Yeckel, Mark Laubach, George K. Aghajanian, and Ralph J. DiLeone (2010) Regulation of Nucleus Accumbens Activity by the Hypothalamic Neuropeptide Melanin-Concentrating Hormone. *J. Neurosci.*, 30: 8263 - 8273.
- Heike Rebholz, Akinori Nishi, Sabine Liebscher, Angus C. Nairn, Marc Flajolet, and Paul Greengard (2009) CK2 negatively regulates G $\alpha_s$  signaling. *PNAS*, 106: 14096 - 14101.
- Emanuela Santini, Myriam Heiman, Paul Greengard, Emmanuel Valjent, and Gilberto Fisone (2009) Inhibition of mTOR Signaling in Parkinson's Disease Prevents L-DOPA-Induced Dyskinesia *Sci. Signal.*, Jul 2009; 2: ra36.
- Zhaoqing Zheng and Joyce Keifer (2009) PKA Has a Critical Role in Synaptic Delivery of GluR1- and GluR4-Containing AMPARs During Initial Stages of Acquisition of In Vitro Classical Conditioning. *J Neurophysiol*, 101: 2539 - 2549.
- Douglas A. Meyer, Edmond Richer, Stanley A. Benkovic, Kanehiro Hayashi, Janice W. Kansy, Carly F. Hale, Lily Y. Moy, Yong Kim, James P. O'Callaghan, Li-Huei Tsai, Paul Greengard, Angus C. Nairn, Christopher W. Cowan, Diane B. Miller, Pietro Antich, and James A. Bibb (2008) Striatal dysregulation of Cdk5 alters locomotor responses to cocaine, motor learning, and dendritic morphology. *PNAS*, 105: 18561 - 18566.
- Akinori Nishi, Mahomi Kuroiwa, Diane B. Miller, James P. O'Callaghan, Helen S. Bateup, Takahide Shuto, Naoki Sotogaku, Takaichi Fukuda, Nathaniel Heintz, Paul Greengard, and Gretchen L. Snyder (2008) Distinct Roles of PDE4 and PDE10A in the Regulation of cAMP/PKA Signaling in the Striatum. *J. Neurosci.*, 28: 10460 – 10471.
- Kurtis D. Davies, Susan M. Goebel-Goody, Steven J. Coultrap, and Michael D. Browning (2008) Long Term Synaptic Depression That Is Associated with GluR1 Dephosphorylation but Not  $\alpha$ -Amino-3-hydroxy-5-methyl-4-isoxazolepropionic Acid (AMPA) Receptor Internalization. *J. Biol. Chem.*, 283: 33138 - 33146.
- Sun X, Milovanovic M, Zhao Y, Wolf ME. (2008) Acute and chronic dopamine receptor stimulation modulates AMPA receptor trafficking in nucleus accumbens neurons cocultured with prefrontal cortex neurons. *J Neurosci*. Apr 16;28(16):4216-30.
- Hu, H., Real, E., Takamiya, K., Kang, M.G., Ledoux, J., Huganir, R.L. and Malinow, R., 2007. Emotion enhances learning via norepinephrine regulation of AMPA-receptor trafficking. *Cell*, 131(1), pp.160-173.

## General References:

- Banke TG, Bowie D, Lee HK, Huganir RL, Schousboe A, Traynelis SF (2000) Control of GluR1 AMPA receptor function by cAMP-dependent protein kinase. *J Neurosci* 20:89-102.
- Esteban JA, Shi SH, Wilson C, Nuriya M, Huganir RL, Malinow R (2003) PKA phosphorylation of AMPA receptor subunits controls synaptic trafficking underlying plasticity. *Nature Neurosci* 6:136-143.
- Hollmann M, Heinemann S (1994) Cloned glutamate receptors. *Annu Rev Neurosci* 17:31-108.
- Keinänen K, Wisden W, Sommer B, Werner P, Herb A, Verdoorn TA, Sakmann B, Seeburg PH (1990) A family of AMPA-selective glutamate receptors. *Science* 249:556-560.
- Lee HK, Takamiya K, Han JS, Man HY, Kim CH, Rumbaugh G, Yu S, Ding L, He C, Petralia RS, Wenthold RJ, Gallagher M, Huganir RL (2003) Phosphorylation of the AMPA receptor GluR1 subunit is required for synaptic plasticity and retention of spatial memory. *Cell* 112(5):631-643.
- Roche KW, O'Brien RJ, Mammen AL, Bernhardt J, Huganir RL (1996) Characterization of multiple phosphorylation sites on the AMPA receptor GluR1 subunit. *Neuron* 16:1179-1188.
- Xiu Sun, Michael Milovanovic, Yun Zhao, and Marina E. Wolf (2008) Acute and Chronic Dopamine Receptor Stimulation Modulates AMPA Receptor Trafficking in Nucleus Accumbens Neurons Cocultured with Prefrontal Cortex Neurons. *J. Neurosci.*, 28: 4216 - 4230.